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Operator's Manual
Manuel d'utilisation
Bedienungsanleitung
Manual del operador

ARGOS
TECHNOLOGIES *manufacturing solutions for science*

Argos Technologies, Inc.
1551 South Scottsdale Court
Suite 200
Elgin, IL 60123
Toll Free: 800-886-8675
International: 847-622-0456
www.argos-tech.com

Table of Contents

Safety Precautions	2
Introduction	3
Product Features	3
Package Contents	3
Specifications	4
Operating Instructions	5
Changing Rotors	6
Cleaning	6
Troubleshooting	7
Warranty	7
Contact Information	7

Safety Precautions

- Do not operate the centrifuge without a rotor properly attached to the shaft.
- Do not fill tubes while they are in the rotor. Liquid spillage may harm unit.
- Do not attempt to remove samples unless the rotor is completely stopped.
- Do not move the centrifuge while the rotor is spinning.
- Do not use a power source other than which the unit designed to operate on.
- Do not use multi-point extension cords with this centrifuge.
- Do not operate centrifuge near a heat source or at temperatures below 4° C.
- Do not open motor casing of centrifuge. This unit is not designed to be user serviceable, and opening by unauthorized personnel will void all warranties.
- Never use solvents or flammables near this or any other electrical equipment.
- Never centrifuge flammable, explosive, or corrosive materials.
- Never centrifuge hazardous materials outside of a hood or proper containment facility.
- Never operate the centrifuge with only one strip tube in place.
- Never operate the centrifuge with uneven loads as this may cause abnormal wear and tear to the motor.
- Always load the rotor symmetrically. Each tube should be counterbalanced by another tube. When spinning 1, 3, 5, or 7 tubes, load a water-filled tube in the opposite slot. When using 0.2mL PCR tubes, load the tubes in diagonally opposite positions to balance the rotor.
- Always install centrifuge in a well ventilated location.
- Always operate centrifuge on a flat surface.
- Always use Allen wrench (provided) to remove the rotor.
- Always locate the centrifuge within easy access to an electrical outlet.
- Always unplug from power before cleaning. Do not immerse.
- Always use proper disinfection procedures when centrifuging biohazardous compounds.
- CAUTION: This unit is not explosion proof.
- CAUTION: Minimize exposure to UV light.

Introduction

The **Argos Technologies Flexifuge Mini Centrifuge** is a benchtop centrifuge that is ideal for quick spin downs from the walls of tubes, microfiltration of HPLC samples, and cell separations. The Flexifuge comes complete with a variety of rotors and features two operating speeds; timed, continuous, and quick-spin modes; and colorful LED lights. The Flexifuge is currently the only microcentrifuge on the market that accommodates 5mL microtubes. With multiple large centrifuge features packed into a personal benchtop unit, the Flexifuge is a cost-effective solution for laboratories on a budget.

Product Features

- Small footprint takes up minimum bench space
- The only microcentrifuge on the market that accommodates 5mL microtubes
- Rotors are easily interchanged by user
- Centrifuge automatically shuts off if top is opened to ensure user safety
- May be used in timed, continuous, or quick-spin modes
- Colorful LED lights alert user at a glance when centrifuging is complete
- Nonskid rubber feet on base prevent vibration and reduce noise levels
- Clear polycarbonate lid allows user to safely observe samples at all times
- Industrial ABS construction is durable and easy to clean


Package Contents

Flexifuge Centrifuge	1 Unit
Cat. Number C1000 (U.S.), C1000E (Europe), C1000U (U.K.)	
Rotor for 1.5/2.0mL tubes	1 Unit
Cat. Number C0300	
Adapters for 0.5mL tubes.	8 Units
Cat. Number C0303	
Rotor for 8-Strip 0.2mL PCR Tubes	1 Unit
Cat. Number C0301	
Rotor for 5mL tubes.	1 Unit
Cat. Number C0302	
Allen wrench	2 Units
Power Supply	1 Unit
Operator's Manual	1 Unit

Specifications

Speed	Slow: 6,000 RPM Fast: 9,200 RPM		
Relative Centrifugal Force		6000rpm	9200rpm
	1.5ml Rotor	2012G	4731G
	0.2ml PCR Strip	1509G	3549G
	5ml Rotor	2414G	5677G
Noise Level	<65dB		
Rotor Capacity	Circular: 8 x 1.5mL tubes (8 x 0.5 with supplied adapters) Strip: Two 8 x 0.2mL tube strips or 16 x 0.2mL tubes Microtube: 4 x 5mL microtubes		
Speed Control	Variable		
Power Requirements	100-240V ~1.0A; 50/60Hz (120V plug supplied) 40 Watts, 1.6A		
Spin Modes	Continuous, Timed, and Quick-Spin		
Timer	1 to 15 Minutes		
Material	Lid: Clear polycarbonate Body: Industrial ABS resin Rotors: Polycarbonate Feet: Silicon Rubber		
Dimensions (L x W x H with lid closed)	9.45 x 7.36 x 5.47 in. (240 x 187 x 139mm)		
Weight	3 lb. 6 oz. (1.6Kg)		
Certifications	CE Marked		

Operating Instructions

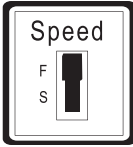1. Select speed by using switch on bottom of unit. Select F for 9,200 RPM or S for 6,000 RPM
2. Select LED color by using switch on bottom of unit: R=Red, B=Blue, G=Green, O=Off
3. Place the centrifuge on a flat, level surface
4. Plug the cord into a properly grounded, three-pronged outlet
5. Open lid by gently pressing the Lid Open Button

2 Tubes


4 Tubes

6 Tubes

8 Tubes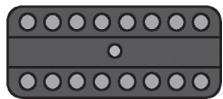

Rotor for 8-strip
PCR Tubes

6. Place tubes into rotor, taking care to load tubes symmetrically

CAUTION: Loading tubes symmetrically ensures quiet operation and significantly extends the life of the centrifuge

7. Set ON/OFF switch to "1" position

Operating Instructions (continued)

8. Make sure that lid is firmly closed. Flexifuge will not operate if lid is not properly latched


9. To use Timed Mode, set Timer to desired length of spin (1 to 15 minutes) and activate by pressing Timer Button

10. To use Continuous Mode, press Continuous Button. To deactivate, press Continuous Button again
11. To use Quick-Spin Mode, press and hold the Pulse Button for desired length of spin. Release Pulse Button to deactivate
12. Wait for rotor to come to a complete stop before removing samples

Changing Rotors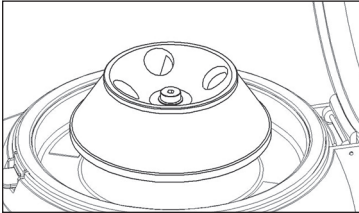

1. Using the Allen wrench loosen the top screw by turning a quarter turn counterclockwise.
2. Remove rotor.
3. Install new rotor and use Allen wrench to tighten locking screw by rotating clockwise.

CAUTION: Operating the 5ml rotor in the fast mode for extended periods of time (>15 minutes) will increase the temperature within the centrifuge.

Cleaning

- Always disconnect unit prior to cleaning.
- To clean the Flexifuge, use a damp cloth and a mild, non-corrosive detergent.
- After cleaning, ensure that all parts are thoroughly dry before attempting to operate the unit.
- Do not immerse the Flexifuge in liquid.

Troubleshooting

Error	Cause	Solution
Unit does not start	Power cable not connected	Check that power cable is firmly plugged into both wall outlet and device
	Power supply	Check that there is power to the wall outlet
	Lid not closed properly	Press firmly to close lid
	Power switch	Check that power switch is at the ON position
Excessive noise or vibration	Rotor is not loaded symmetrically	Stop the unit and load the tubes symmetrically. Ensure that the tubes on the opposite sides are equally filled.
	Rotor is slanted	Remove the rotor and reinsert straight and firmly downwards
	Loose locking screw	Check to see if locking screw has been inserted properly and that it has not been over-tightened
	Unit base is not level	Check that unit on a flat surface
If suggested solutions are unsuccessful, please contact Argos Technologies at 847-622-0456		

Warranty

- The full warranty for this device is 12 months from the date of sale.
- During this period of time the device will be repaired or exchanged without charge.
- Warranty is null and void on devices with damaged manufacturer's number; damaged seal; damage due to improper use; storage, or transport of device; damage due to service by unauthorized persons.

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Copyright © 2009 Argos Technologies, Inc.