Argos Technologies RotoFlex™ Tube Rotator

Operator's Manual

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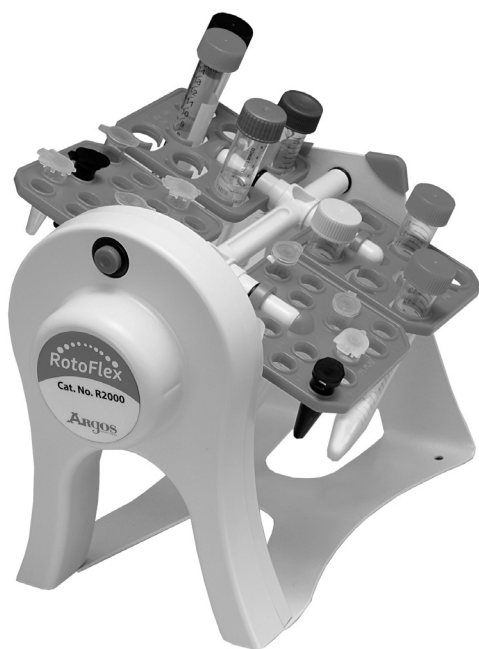
Manuale d'uso

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ARGOS
TECHNOLOGIES *manufacturing solutions for science*

Argos Technologies, Inc.

1551 South Scottsdale Court

Suite 200

Elgin, IL 60123

Toll Free: 800-886-8675

International: 847-622-0456

www.argos-tech.com

Introduction

Thank you for purchasing the **Argos Technologies RotoFlex Tube Rotator**. The unit was designed to provide researchers the ability to mix a variety of tube sizes (.5 - 50mL) with a 360° rotation. The adjustable rotisserie also allows you to set the mixing angle from 0-90°; depending on the application.

To change rotisserie, simply snap on the desired tube size to the support bar. No tools required.

The RotoFlex is ideal for

- Hybridizations
 - DNA purification
 - Sample mixing / resuspensions
 - Protein binding
 - Immunoprecipitation
 - Any procedure where thorough mixing is required
-

Package Contents

Tube Revolver.	1 Unit
Cat. Number R2000 (120V), R2001 (220V)	
Rotisserie for 0.5 - 1mL tubes	1 Unit
Cat. Number R2002	
Rotisserie for 1.5 - 2mL tubes	1 Unit
Cat. Number R2003	
Rotisserie for 5 - 7mL tubes.	1 Unit
Cat. Number R2004	
Rotisserie for 10 - 15mL tubes.	1 Unit
Cat. Number R2005	
Rotisserie for 50mL tubes	1 Unit
Cat. Number R2006	
Revolving Rack.	1 Unit
Operator's Manual	1 Unit

Specifications

Model	R2000	R2001
Rotation Speed	22rpm	18rpm
Capacity	42 x 0.5-0.8mL tubes / 30 x 1.5-2.0mL tubes / 14 x 5-7mL tubes / 14 x 10-15mL tubes / 6 x 50mL tubes	
Ambient Temperature	4°C to 60°C 20-85% RH non-condensing	
Speed Control	Fixed	
Max Load	1.8lb / 0.8kg (Comprise Rotisseries)	
W x D x H (cm)	22.5 x 15.4 x 19.5	
Power	110 V 60 Hz	220 V 50Hz
Noise Level	Noisless Operation	
Weight	1.8 lb / 0.8kg	
Certifications	CE Marked	

This device complies with RoHS 2009/95/EC

Operating Instructions

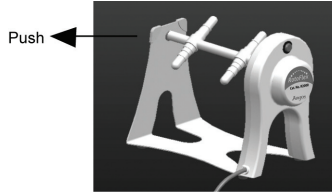
1. Make sure the power switch is turned off before connecting the plug to a grounded wall socket.
2. Place the RotoFlex on a level, horizontal and secure surface.
3. Select the rotisserie required for your application (see page 4) for rotisserie installation.
4. Load the tubes and adjust the rotisserie to the desired angle. Be sure to distribute tubes evenly to achieve a smooth rotating motion.
5. Start the rotator by pushing down the power switch.
6. When finished, turn the power switch off before removing tubes.

Note: You can change the rotation direction by pushing the rotisserie in the opposite direction.


Caution: Do not touch plug with wet hands or use the device near a water source.

Changing Rotisseries

1. Firmly grasp the RotoFlex rotator with your left hand while holding the rotisserie axle with your right hand. Gently pull the rotator body with your left hand outward until the axle is free of the unit.

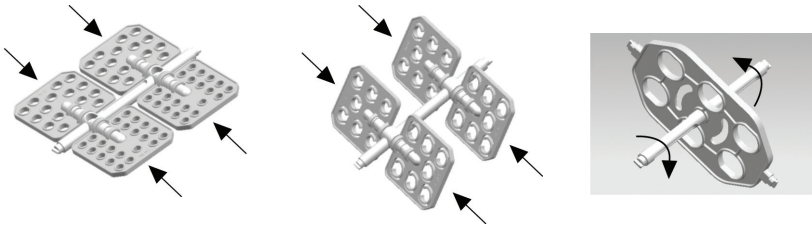2. Remove the opposite side of the axle body by pulling it gently from the RotoFlex housing.
3. Select the rotisserie you wish to utilize and mount them on the RotoFlex axle.

4. Insert the axle back into the RotoFlex housing and turn until the axle clicks into place.
5. Gently pull the rotator body (as in step 1) and reinsert the axle.
6. Turn the unit on to allow the rotisserie axle to align with the RotoFlex motor.

Maintenance & Cleaning

- No lubrication is required.
- Before cleaning be sure to unplug the power cord from the wall socket.
- Remove the rotisseries.
- The housing can be cleaned using a soft damp cloth and a mild non-corrosive detergent.
- After cleaning, ensure that all parts are thoroughly dry before attempting to operate the unit.
- **Do not immerse the RotoFlex in liquid.**

Warnings

- Not to be used to mix flammable samples.
- Avoid aggressive solvents or vapors.
- Take appropriate precautions when mixing hazardous samples.

Troubleshooting

Error	Cause	Solution
Unit does not start	Power cable not connected	Check that power cable is firmly plugged into wall outlet.
	Power Supply	Check that there is power to the wall outlet.
	Power Switch	Check that power switch is at ON status.
<p>If suggested solutions are unsuccessful, please contact Argos Technologies Technical Support at</p> <p>Argos Technologies 800-886-8675 (Inside U.S.) 847-622-0454 (Outside U.S.) www.argos-tech.com</p>		

Warranty

Argos RotoFlex Tube Rotator is warranted to be free of defect in materials and workmanship for a period of 12 months from the date of purchase, providing that:

- The unit has not been opened or attempted to be opened by the end user.
- The user ensures that it has been carefully cleaned and decontaminated.
- The unit has been operated according to the instructions and not in any improper way.

In order to qualify for warranty repair; the customer must return the warranty card supplied with the unit that includes all of the information required to validate proof of purchase. Should warranty service be required, repair or replacement will be provided by Argos without charge. Freight shipping the unit to Argos will be the responsibility of the purchaser. There are no other warranties or guarantees, implied or otherwise expressed including any warranties of application or merchant

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Copyright © 2011 Argos Technologies, Inc.